

H 318 Causes serious eye damage.
P 280-305+351 +338-310

Nitrite No. 1



H 411 Toxic to aquatic life with long lasting effects.
P 273-391-501

Nitrite No. 2



Water-i.d./Pool-i.d. GmbH
Daimlerstr. 20
D-76344 Eggenstein
Tel. +49 (0) 721 - 782029-0
Fax +49 (0) 721 - 782029-11
www.water-id.com
info@water-id.com

NOTRUF/EMERGENCY: +49 (0)89-19240 (24h)



Nitrite No. 1

CORROSIVE



Nitrite No. 2

**DANGEROUS
FOR THE ENVIRONMENT**



Nitrite

Nitrit

Nitrite

Nitrito

Nitrito

Nitrito

Nitrit

Nitriet

Tablet-count method

Tabletten-Zählverfahren

Methode de comptage à pilule

Metodo di conteggio de pastiglie

Método Contador de tabletas

Método de contagem de comprimidos

Tablettællings-metoden

Tablet-tælmethode

Range: 70 - 700 mg/l NaNO_2 : 20 ml sample water + 30 ml aqua dest!


Range: 140 - 1400 mg/l NaNO_2 : 10 ml sample water + 40 ml aqua dest!

Range: 280 - 2800 mg/l NaNO_2 : 5 ml sample water + 45 ml aqua dest!




①

Fill in
sample
water!

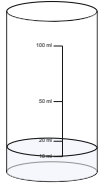


②

Add
aqua
distilled!

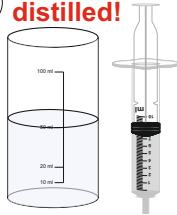


③

Nitrite No. 1

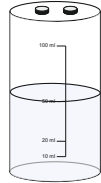


④

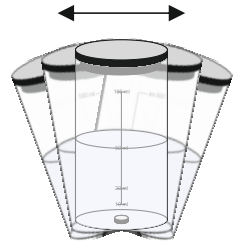


⑤



Nitrite No. 2

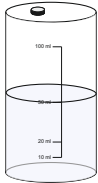

⑥

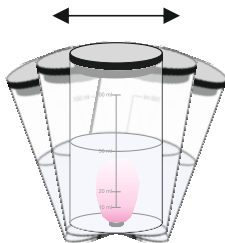

⑦

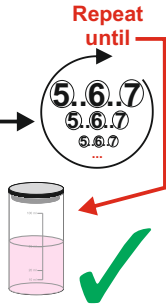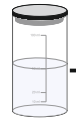

Range 70 - 700 mg/l:  Nitrite No. 2 x 70 =

Range 140 - 1400 mg/l:  Nitrite No. 2 x 140 =

Range 280 - 2800 mg/l:  Nitrite No. 2 x 280 =



QUALITY REAGENTS



MADE IN GERMANY